



# PROMED

A PARTNERSHIP IN PRACTICE

## Lec Control Plus

### Checking and Resetting Your Recorded Min/Max Air Temperatures



#### STEP 1

##### ENSURE AIR TEMPERATURE IS SELECTED

The Air/Load button toggles between the air and load temperatures.

Please ensure "A" is displayed on the top right of the screen before checking and recording temperatures.

If "A" is not displayed, press the Air/Load button

Further information on the Load temperature can be found in the fridge manual



#### STEP 2

##### View the maximum recorded temperature

Press the MAX button

The screen will show:

- The word HI for 2 seconds
- Then the **maximum recorded** temp. for 2 seconds
- Then returns to the running air temp.



#### STEP 3

##### View the minimum recorded temperature

Press MIN button

The controller shows

- The word LO for 2 seconds
- Then the **minimum recorded** temp. for 2 seconds
- Then returns to the running air temp.



#### STEP 4

##### To RESET BOTH the AIR and LOAD

MIN and MAX temperatures

Press and hold the RESET button for 5 seconds

Once the reset is complete, the controller will beep

Call us on Freephone (ROI) 1800 619 619

# LEC Fridge

## Correct stocking example

For optimum airflow, do NOT fill your refrigerator above 75% capacity, and leave gaps between products. Ensure the door is opened for minimal time when restocking

Products **NOT** blocking the fan for optimal airflow



Products not stacked against the back or side walls

Space between products to aid airflow



Products stored in the basket and not on the base of the fridge